



# SWIFT

OPTICAL INSTRUMENTS

Exceptional Optics ■ Durable Construction ■ Innovative Designs

## Wired Digital Microscope or Camera Connected to Laptop Connected Wirelessly to Router



Start with a wired digital microscope or camera connected to Motic Images Plus software. MoticHUB is built into the software. The teacher's WiFi enabled laptop must be wirelessly connected to the router.

Students using tablets or smart phones can connect to the digital microscope or camera using the free MotiConnect app and the IP address generated by MoticHUB.



Students connected to the network and using Motic Images Plus (must be installed on the local drive) can connect to the digital device by inputting the IP address generated by MoticHUB.